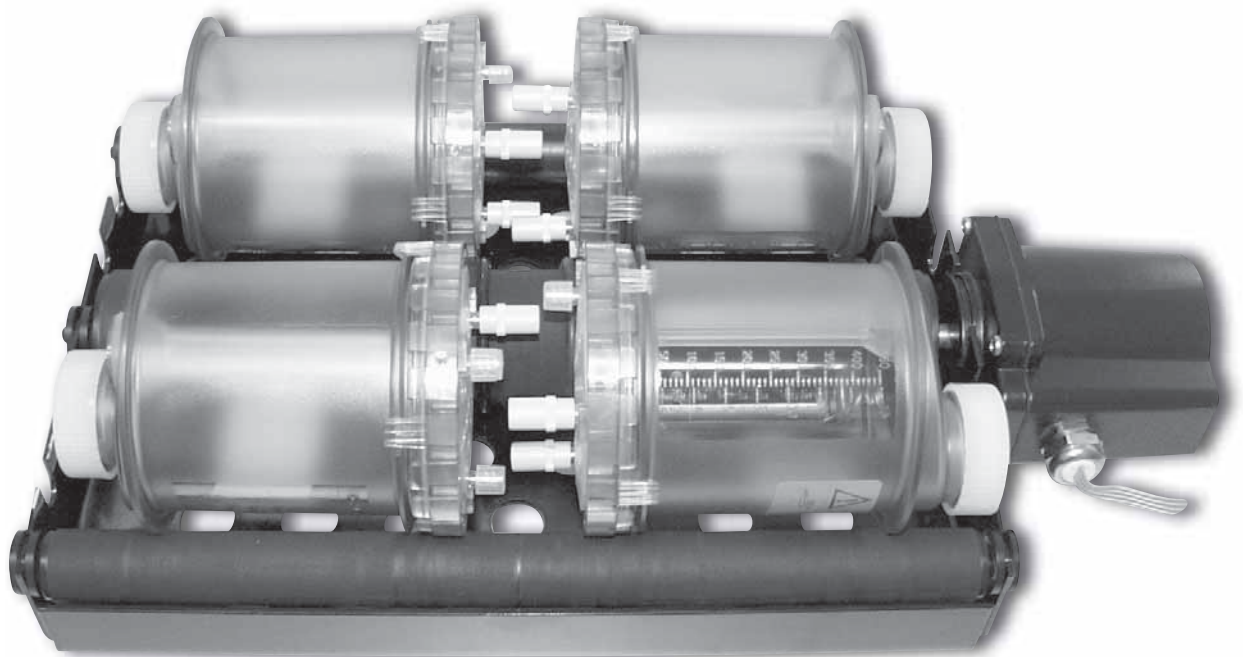
---

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Universaldrehvorrichtung

CE

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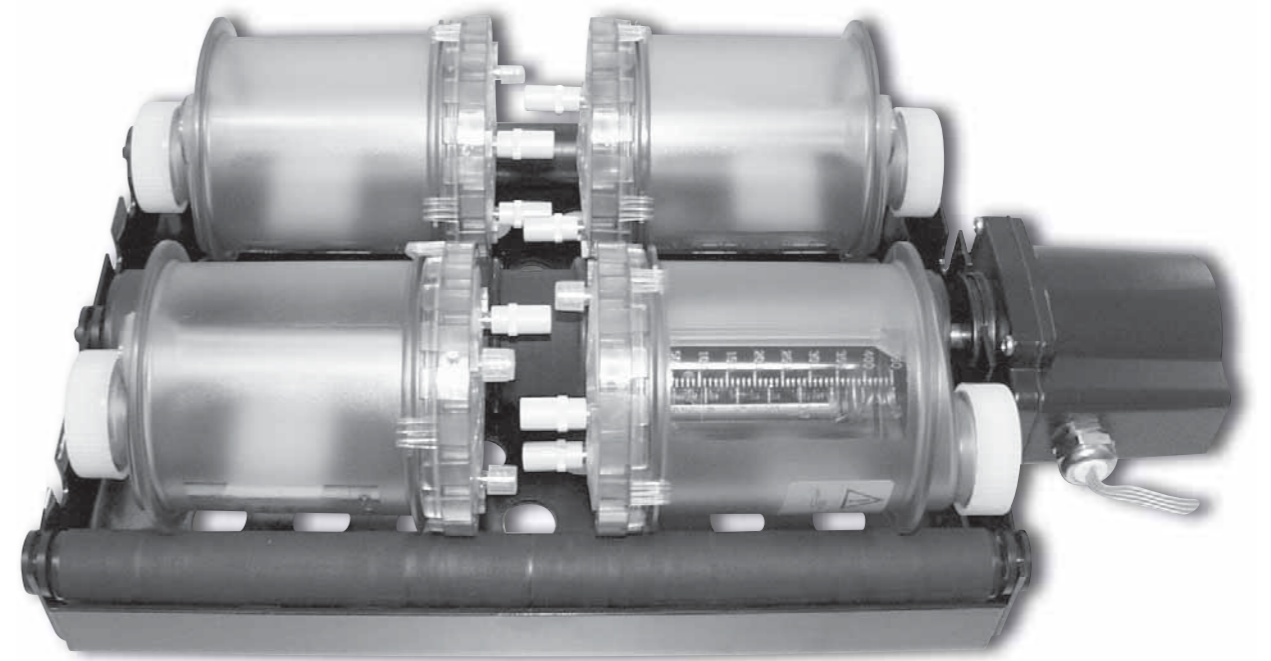
3. Important instructions ..... 3

    Explanation of symbols ..... 4

    Device-specific safety instructions..... 4

    Operator obligations ..... 5

    Staff obligations ..... 5

4. Further instructions ..... 5

    Guarantee and liability ..... 5

    Decommissioning/ Disposal ..... 6

5. Description of the device..... 6

6. Transport, installation and connection ..... 6

    Scope of delivery ..... 6

    Unpacking ..... 7

    Specifications ..... 7

    Installation ..... 7

    Startup ..... 7

7. Operation ..... 8

    Controls ..... 8

8. Service and maintenance ..... 10

    Cleaning ..... 10

    Service address ..... 10

    Contamination questionnaire..... 11

**1. Device data**

*(to be completed by the customer)*

Type: \_\_\_\_\_

Serial No.: \_\_\_\_\_

Place of installation: \_\_\_\_\_

Installation date: \_\_\_\_\_

Inventory No.: \_\_\_\_\_

**2. Manufacturer's address**

SARSTEDT AG & Co.  
P.O. Box 12 20  
D-51582 Nümbrecht  
Phone +49 2293 3050  
Fax +49 2293 305- 122  
info@sarstedt.com

**3. Important instructions**

**Follow this user guide's safety instructions before you first operate the device!**

Awareness of the content of this User Manual is a prerequisite for correctly handling the device and for smooth operation.

Please keep the User Manual to your device for future consultation.

**Copyright:**

SARSTEDT AG & Co owns the copyright to this User Manual. The User Manual is only intended for use by the staff operating the device and the purchaser. This User Manual must neither be reproduced in full or in part nor distributed without the prior written consent of SARSTEDT AG & Co. Any infringements will be prosecuted.

Technical modifications reserved!

Nümbrecht, November 2009

SARSTEDT AG & Co.

### Explanation of symbols

This User Manual includes the following important symbols to indicate hazards and operating errors.

The „warning“ symbol means that this section:

- contains important instructions.
- describes an immediate or potentially lethal danger or threat to a person's health.
- contains a reference to dangerous situations.

**Ignoring this symbol can cause light to serious or lethal injuries or damage to property.**

**Warning of dangerous electrical voltage!**

This „information“ symbol refers to information on the proper handling of the device or to general information to facilitate understanding. Failure to observe these instructions may lead to improper operation or damage to the device.

Other symbols refer to the topic depicted in the symbol, such as to regulations for the prevention of accidents, power supply, servicing or disposal.

### Device-specific safety instructions

The device has been designed in accordance with state-of-the-art technology and recognised safety-specific regulations. The appropriate safety instructions and guidelines in addition to the occupational safety and accident prevention procedures for use in laboratories must be duly observed.

Nevertheless, danger to the user, third-parties or damage to the device or other property can occur during operation.

Only use the device:

- for its intended purpose.
- in perfect condition with regard to safety.

The device must not be operated in explosive areas or used for rolling explosive or flammable substances. The device is a laboratory instrument and conforms to the international laboratory device safety regulations (IEC 61010).

### Device-specific safety instructions

The device can be operated at a voltage of 115 or 230 V/50-60 Hz.

It is, therefore, vital to adhere to the following instructions:

- Do not let any water or other liquids enter the device.
- Only use the adaptor plug provided to connect the device to the mains.
- Check the connection cable for damage before operation.
- Never use any defective cables.
- Only connect the device to the mains voltage indicated on the rating label.
- Do not operate defective devices or cables.
- Never open parts of the casing held together by screws.
- Unplug the adaptor from the socket for cleaning or if the device is not in use for any length of time.

Only attach accessories or operate the device with those specifically suited for these purposes.

### Operator obligations

The operator undertakes to assign work on the device only to those individuals who have read and understood this User Manual. This should be confirmed by signature. Staff awareness as to safe operation should be checked at regular intervals.

### Staff obligations

Before initial startup, individuals working with the device for the first time undertake to:

- read this User Manual.
- follow the prevailing rules and procedures for accident prevention.

### 4. Further instructions

#### Guarantee and liability

As a rule, the 'General conditions of sale' of SARSTEDT AG & Co. apply. They are listed on the back of the invoice. During the guarantee period the device may only be repaired by SARSTEDT AG & Co. or by individuals authorized by SARSTEDT AG & Co.. Any warranty claim will expire in case of improper handling or repair. Warranty and liability claims are excluded if operation of the device inconsistent with its intended use caused by one or more of the following reasons:

- operation of the device inconsistent with its intended use.
- improper assembly, initial startup, operation and maintenance of the device.
- device operation with defective safety equipment or improperly attached or non-functioning safety and protective devices.
- non-observance of the instructions in the User Manual on transport, storage, assembly, startup, operation, servicing, setting-up and disposal.
- unauthorised modifications to the device.
- emergencies resulting from contact with foreign bodies and force majeure.
- improperly performed repairs.

### Decommissioning / Disposal

- It is imperative to handle the device professionally and properly and dispose of it in accordance with the effective legal regulations.
- It is imperative to handle any substances used in connection with this device professionally and properly and to dispose of them in accordance with the effective legal regulations.

This user guide consists of the following materials:  
The back of the binder and sheet cover are made of PVC, the remainder is made of paper.  
The paper can be separated from the plastic waste by pulling off the back of the binder.

### 5. Description of the device

The Universal Turning Device is used for cultivating cells in the miniPERM® bioreactor. Designed for up to four bioreactors (max. loading weight 3 kg) and an individually adjustable rotational speed (0.1 - 40 rpm), the device provides optimum requirements for each cell type. The control unit for the Universal Turning Device is placed and also operated outside the incubator, and therefore particularly space-saving. An extensive range of accessories enables the device to be used multi-functionally in addition to cell cultivation. In order to expand the possible applications, we provide adaptors for centrifuge tubes and reagent vessels and also a rocker device.

### 6. Transport, installation and connection

#### Scope of delivery

##### Standard delivery includes

1x Tray  
2x Side rollers  
1x Drive roller  
4x Spacers  
(on the drive rollers)  
1x Motor  
1x Control unit  
2x Fuses (160mA, slow-acting)

##### Accessories (optional)

Art. No.: 94.6077.360  
Adaptor for centrifuge tubes  
12 x 15 ml  
Art. No.: 94.6077.361  
Adaptor for centrifuge tubes  
5 x 50 ml  
Art. No.: 94.6077.362  
Adaptor for reagent vessels  
48 x 1.5/2.0 ml  
Art. No.: 94.6001.000  
Rocking device top part set

##### Spare parts

Art. No.: 94.6291.061  
Control unit  
Art. No.: 94.6001.064  
Motor  
Art. No.: 94.6055.594  
Side rollers  
Art. No.: 94.6055.593  
Drive roller  
Art. No.: 94.6055.440  
Tray  
Art. No.: 94.6290.560  
Spacer

If you have any queries relating to SARSTEDT's Universal Turning Device, accessories or the SARSTEDT product range, please contact:

**SARSTEDT AG & Co.**  
P.O. Box 12 20  
D-51582 Nümbrecht  
Phone +49 2293 3050  
Fax +49 2293 305-122  
info@sarstedt.com

or your local SARSTEDT sales organisation.

### Unpacking

- Check the integrity of the cardboard packaging and the device.
- Contact the carrier without delay in the event of transport damage! Observe the notification period of the individual transport companies (railway, post office, parcel service or carrier) which may be only 24 hours.
- Check items delivered
- Report defects and damage immediately to SARSTEDT AG & Co.!
- To prevent damage to individual parts, do not open the packaging with a knife or sharp item.

### Specifications

#### Device data

Device: Universal Turning Device  
Part number: SARSTEDT No. 94.6001.061

#### Power supply

Electrical connection: 115/230 V; 50/60 Hz  
Power consumption: 18 VA

#### Dimensions

Tray (including motor on drive shaft): W x H 46 x 26 cm  
Control unit: W x H 22 x 10 cm

#### Weight

Tray (including rollers and motor): 4.2 kg  
Control unit: 1.4 kg

#### Ambient conditions

Admissible ambient temperature: +15 - +40 °C (motor max. +60 °C)  
Max. relative humidity: 80%

### Installation

- Place the device on a firm vibration-free and level surface.
- Maintain at least 10 cm clearance to other items and devices.
- There must not be any contact with other devices or items during operation.
- The device's power switch must be easily accessible.

### Startup

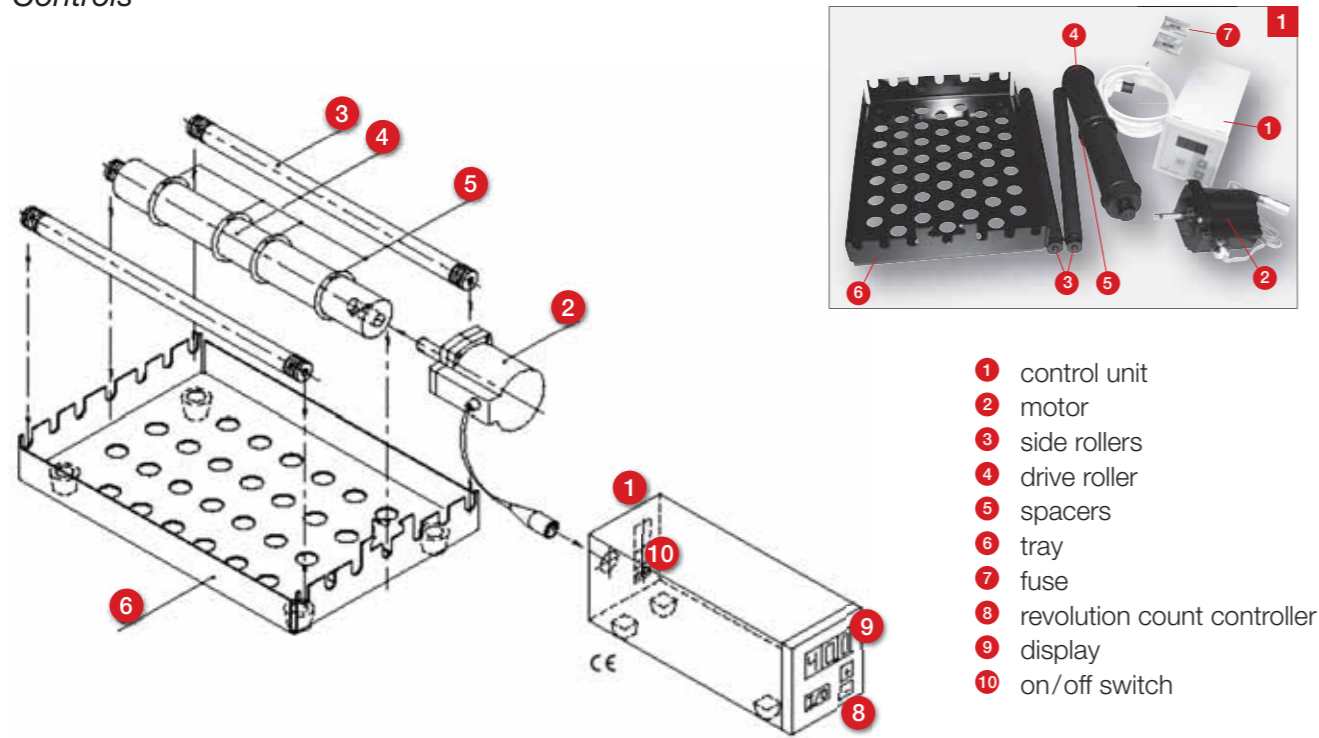
Warning before initial startup:

- Check if the current mains voltage matches the rated voltage (See Safety Instructions, page 9).
- Adhere to the admissible ambient conditions for the device (See Specifications, page 7).
- In the cold season wait approx. ½ to ¾ of an hour after delivery and unpacking for the device to reach ambient temperature (18 °C to 23 °C).


After connecting to the mains, the device is ready for use.

7. Operation  
Controls


- 1 control unit
- 2 motor
- 3 side rollers
- 4 drive roller
- 5 spacers
- 6 tray
- 7 fuse
- 8 revolution count controller
- 9 display
- 10 on/off switch

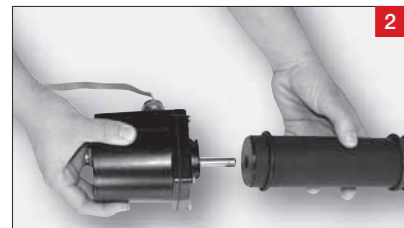Insert the motor drive shaft into the mounting hole of the drive roller.


Note!

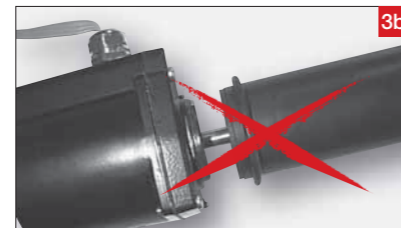The notch on the drive axle must lock into the driving pin of the drive roller (3a/3b).

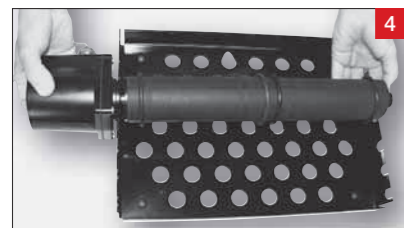
Mount the motor and drive roller into the tray. First mount the drive roller bearing into the mounting device on the tray. Then push the notch on the motor case into the semicircular tray holder and mount it.

Note!

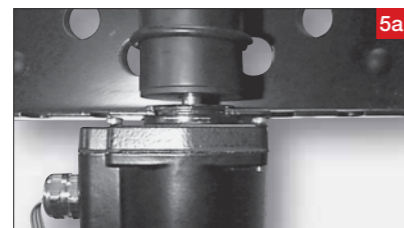The motor case notch must lock into the semicircular tray holder. Do not let the drive axle rest upon the tray holder (5a/5b).

Mount the side roller bearings into the tray holder. (Select the outer openings when using the Universal Turning Device for miniPERM® bioreactors.)

Connect the flat cable on the motor to the control unit. Place the assembled tray into the incubator and the control unit outside the incubator. Run the flat cable through the door seal.

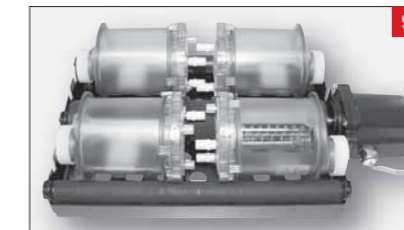
Position the spacers on the drive roller at a sufficient distance to the ends of the roller (min. 30 mm).

If the distance to the ends of the roller, is too small, there is the risk of the miniPERM® bioreactors jamming on the tray during rotation (8a/8b).

When using miniPERM® bioreactors with in-stoppers instead of cylindrical screw caps, the bioreactors must be positioned on the universal turning device as in Fig. 9. The two central drive roller spacers require at least 25 mm clearance.

Safety Note

**Do not connect the mains cable beforehand!**

The Universal Turning Device, as-delivered, is preset to a voltage of 230 V (control window: 230 V, fig. 10). Observe the following instructions in case of deviating voltage (115 V):

10 Adjust the voltage switch to the required voltage (factory setting: 230 V, 2 x 80 mA, slow-acting).

11 Open the fuse cover and take the fuse holders out of the device.

12 Remove both fuses and replace with the ones supplied (2 x 160 mA, slow-acting).

13 Turn the fuse holders through 180° (control window: 115 V) and re-insert them.

8. Service and maintenance

Cleaning

**It is imperative to observe the instructions for cleaning the device. Negligent cleaning or ignoring the instructions can cause malfunctions!**

- Disconnect the device to be cleaned from the power supply.
- Do not let any liquid enter the device while cleaning.
- Clean parts of the casing with a slightly damp cloth. You can use 70% ethanol. Dry the device carefully afterwards.
- Never use scouring agents, caustic cleaners or solvents (except alcohol).

**You can clean and disinfect each part of the Universal Turning Device with 70% ethanol.**

Service address

If you have any queries or problems concerning the device, please contact the DESAGA GmbH/SARSTEDT GROUP service department in Wiesloch or your local SARSTEDT sales organisation. Always state the device's serial number.

**DESAGA GmbH / SARSTEDT GROUP**  
**Serviceabteilung**  
**In den Ziegelwiesen 1-7**  
**D-69168 Wiesloch**  
**Phone: +49 62 22 / 92 88 65**  
**Fax.: +49 62 22 / 92 88 60**

- Complete the contamination questionnaire (see the end of the user guide) and include with the device a copy of the delivery note and brief details of the problems that have occurred.
- Always ship the device in its original packaging. If it is no longer available, use packaging which will prevent damage to the device during transport. The sender assumes responsibility.
- The contamination questionnaire is for the safety of our service staff.

Please complete it as carefully as possible in order to avoid delays in repair processing.

**Contamination Questionnaire for Repair Work**

**Dear customer,**

For safety reasons please answer all the following questions before dispatching your device for service, repair or return to DESAGA GMBH/SARSTEDT GROUP, Wiesloch.

Company: \_\_\_\_\_ Department: \_\_\_\_\_

Town: \_\_\_\_\_ Street: \_\_\_\_\_

Name: \_\_\_\_\_ Telephone No.: \_\_\_\_\_

Device/Art. No.: \_\_\_\_\_ SN: \_\_\_\_\_

- The device is free of hazardous materials
- The device was in contact with the following hazardous materials

	Material class	Material name
<input type="checkbox"/>	Corrosive materials	
<input type="checkbox"/>	Easily inflammable materials	
<input type="checkbox"/>	Environmentally hazardous materials	
<input type="checkbox"/>	Poisonous materials	
<input type="checkbox"/>	Harmful materials	
<input type="checkbox"/>	Radioactive substances	
<input type="checkbox"/>	Biologically hazardous materials	
<input type="checkbox"/>	Other hazardous materials	
<input type="checkbox"/>	The device has been decontaminated according to legal regulations	
(Tick where applicable)		

**Description of decontamination work performed:**

Date: _____	Signature: _____

Technical modifications reserved

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