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SARSTEDT Transferpipetten

DE	Gebrauchshinweis – Transferpipetten	2 - 3
EN	Instructions for Use – Transfer pipettes	4 - 5
BG	Инструкции за употреба – Пастър пипети	6 - 7
CS	Návod k obsluze – Transfer pipety	8 - 9
DA	Brugsanvisning – Transfer pipette	10 - 11
EL	Οδηγίες χρήσεως – Πιπέτες μεταφοράς	12 - 13
ES	Instrucciones de uso – Pipetas de transferencia	14 - 15
ET	Kasutusjuhend – Ülekandepipett	16 - 17
FI	Käyttöohje – Pasteur pipetti	18 - 19
FR	Mode d'emploi – Pipettes de transfert	20 - 21
HR	Uputa za upotrebu – Transfer pipete	22 - 23
HU	Használati utasítás – Transzferpipetták	24 - 25
IT	Istruzioni d'uso – Pipette di trasferimento	26 - 27
KO	사용 설명서 – 이송 피펫	28 - 29
LT	Naudojimo instrukcijos – Perkėlimo pipetės	30 - 31
LV	Lietošanas norādes – Pastēra pipetes	32 - 33
NL	Gebruiksaanwijzing – Transferpipet	34 - 35
NO	Bruksanvisning – Transferpipette	36 - 37
PL	Instrukcja obsługi – Pipety transferowe	38 - 39
PT	Instruções de utilização – Pipetas de transferência	40 - 41
RO	Instrucțiuni de utilizare – Pipete de transfer	42 - 43
RU	Инструкция по применению – пипетки Пастера	44 - 45
SK	Návod na Použitie – Transfer pipety	46 - 47
SL	Navodila za uporabo – Pipete za prenos	48 - 49
SV	Bruksanvisning – Transferpipetter	50 - 51
TH	คำแนะนำในการใช้งาน – ปิเปตต์แบบปริมาตร	52 - 53
TR	Kullanım için talimatlar – Transfer pipetleri	54 - 55
ZH	使用说明 – 移液器	56 - 57

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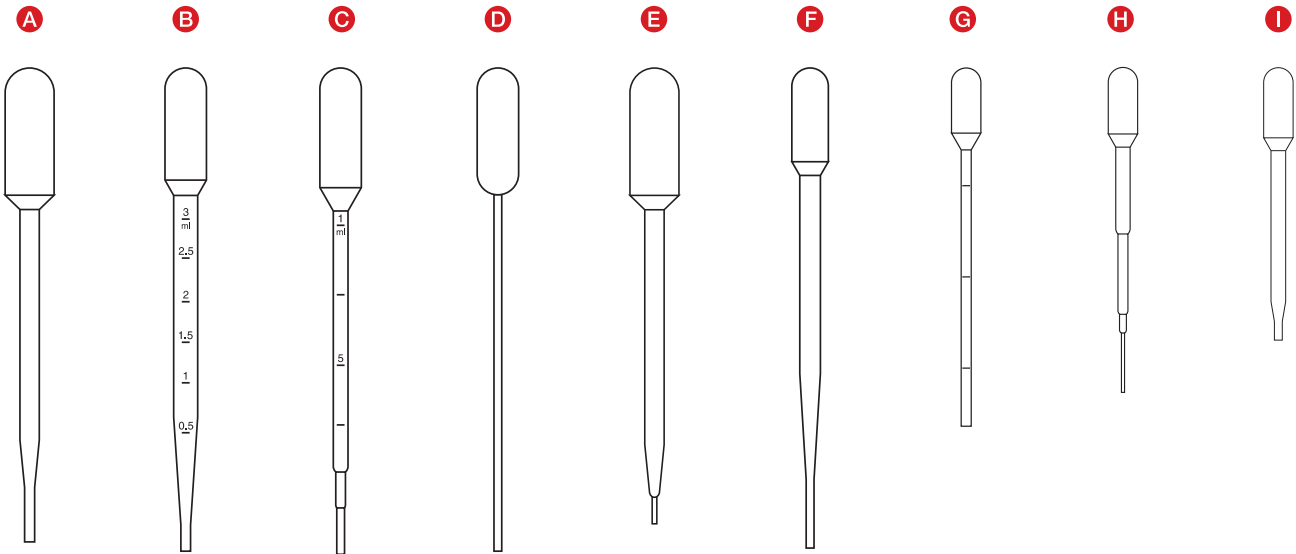
Intended use

Transfer pipettes are used for transferring diagnostic sample material from one vessel to another. This allows fluids such as urine, blood, serum and plasma to be transferred for subsequent analysis. Transfer pipettes are used when it is not necessary for the sample volume to be exact. If exact volumes are needed, piston pipettes with suitable pipette tips should be used. No special ambient conditions are required. Transfer pipettes should be used by medical professionals.

Product description

Transfer pipettes with an integrated suction bulb are available in nine different versions. The different versions (A-I) vary in terms of suction volume, droplet size, length and graduation. In addition, sterile and non-sterile variants are available, as well as different packaging options. Sterile transfer pipettes are used, for example, for transferring sample material for microbial analyses. Here, there is a choice between individual sterile packs and sterile packs of 10 transfer pipettes.

Transfer pipettes:

Type	Art. no.	Name
A	86.1174.XXX	6 ml LD-PE transfer pipette, without graduation, Length: 152 mm, droplet size: 40 - 55 µl
B	86.1171.XXX	3.5 ml LD-PE transfer pipette, with graduation 3: 0.50 ml (0.5 - 3.0 ml), Length: 155 mm, droplet size: 35 - 55 µl
C	86.1172.XXX	3.5 ml LD-PE transfer pipette, with graduation 1: 0.25 ml (0.25 - 1.0 ml), Length: 156 mm, droplet size: 30 - 45 µl
D	86.1173.XXX	3.5 ml LD-PE transfer pipette, without graduation, Length: 155 mm, droplet size: 30 - 45 µl
E	86.1175.XXX	6 ml LD-PE transfer pipette, without graduation, Length: 146 mm, droplet size: 20 - 30 µl
F	86.1176.XXX	2 ml LD-PE transfer pipette, without graduation, Length: 154 mm, droplet size: 35 - 45 µl
G	86.1170.XXX	1 ml LD-PE transfer pipette, with graduation 1: 0.10 ml (0.1 - 0.3 ml), Length: 115 mm, droplet size: 45 - 55 µl
H	86.1180.XXX	1 ml LD-PE transfer pipette, without graduation, Length: 104 mm, droplet size: 17 - 22 µl
I	86.1179.XXX	1 ml LD-PE transfer pipette, without graduation, Length: 87 mm, droplet size: 30 - 45 µl

LD-PE = low density polyethylene

Storage

Store the product at room temperature.

Handling

1. For blister pack variants: Open the blister at the opening tab marked with corrugations. This tab is located on the side of the suction bulb of the transfer pipette. The transfer pipette can be grasped by the suction bulb and removed from the blister pack.
2. Press in the suction bulb of the transfer pipette.
3. Dip the tip of the transfer pipette into the fluid you intend to transfer.
4. Release the suction bulb, allowing the vacuum created to fill the transfer pipette with fluid.
5. Transfer the tip of the transfer pipette into the target vessel and press the suction bulb again to release the fluid.
6. Dispose of the transfer pipette in accordance with the regulations of your facility.

Restriction

- The accuracy of the droplet size and the graduation is $\pm 15\%$. Transfer pipettes are therefore intended for transferring undefined fluid quantities. For pipetting exact volumes, a piston pipette with a suitable pipette tip should be used.

Disposal

- Wearing disposable gloves and other personal protective equipment reduces the risk of infection. Follow the regulations of your facility.
- General hygiene guidelines and legal regulations for the proper disposal of infectious material should be observed and complied with.
- Contaminated transfer pipettes must be disposed of in appropriate biohazard disposal containers. Follow the regulations of your facility.

Safety information/warnings

- Personal protective equipment must be used when handling diagnostic samples and hazardous materials.
- Do not use transfer pipettes after the expiry date has passed
- The handling of biological samples must take place in accordance with the applicable hygiene and safety regulations.
- Transfer pipettes are intended for single use. After use they should be disposed of according to the applicable regulations. There is a risk of contamination if reused.

Product-specific standards and guidelines in 80/181/EEC as amended/updated:

Units in the measuring system

Key for symbols and labels:

Article number

In vitro diagnostic device

Batch number

CE mark

Manufacturer

Store at room temperature

Sterilisation with ethylene oxide gas

If reused: Risk of contamination

Do not resterilise

Use by

Follow the instructions for use

Country of manufacture

Keep away from sunlight

Store in a dry place

Do not use if package is damaged

Single sterile barrier system

Single sterile barrier system with protective packaging outside

Authorised representative in Switzerland

Technical modifications reserved.

All serious incidents relating to the product must be reported to the manufacturer and the competent national authority.

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